



65 Commerce Way Dedham, MA 02026
 Phone 1-800-796-9199 or 781-326-6709
 Fax 1-781-326-5611 Web www.divbio.com

FORMATTING INSTRUCTIONS FOR LASER CRYO-BABIES® and CRYO-TAGS® IN MICROSOFT WORD®

1. Open Word
2. Click on Mailings > Labels
3. Click on Labels tab
4. Click on Options
5. Under Tray, select Default tray (Automatically Select) or Manual Feed
6. Click on New Label to customize your label
7. Type in a name for the label (e.g. Catalog Number)
8. Insert the Label Details **[shown at bottom of page]** and click OK
9. Select the label you made in the Label Options window and click OK
10. In the Envelopes & Labels window, type your information in the area provided and click on New Document
11. When the cells appear on the screen, select the entire document
12. When the whole document is highlighted, choose an appropriate font and font size* and center the type horizontally using the Center command located in the toolbar

If the text in either the rows or columns do not align, continue with Steps 14 -15 otherwise skip to Step 16.

13. After centering the text horizontally, highlight the entire sheet and align the text vertically by right clicking and selecting Cell Alignment
14. Click the Align Top Center, Left or Right icon (depending on where the text needs to be aligned)
15. To determine if your text will fall within the boundaries of the label, have the Gridlines showing
16. Before printing a sheet of labels, print a test page on blank paper and line it up with a sheet of labels to ensure the text lines up correctly

Notes:

- Pass the laser sheet through the printer only once
- Manually feed the laser sheets one at a time
- Print your labels on a blank sheet of paper and check the positioning on a blank sheet of labels
- Minor adjustments may be required depending on your printer and/or software you are using

* Suggested Font Sizes: 8 pt. for smaller labels or 10-12 pt. for larger labels

Catalog No.	LCRY-1700	LCRY-1200	LCRY-2380	LCRY-1100	LCRY-1258	SIDE-1000
Label Size	1.28" x 0.50"	1.50" x 0.75"	0.94" x 0.50"	1.69" x 0.75"	2.625" x 1.00"	1.50" x 0.25"
Top Margin	0.24"	0.31"	0.24"	0.63"	0.50"	0.62"
Side Margin	0.77"	0.25"	0.56"	0.70"	0.19"	0.50"
Vert. Pitch	0.63"	0.88"	0.63"	0.75"	1.00"	0.25"
Horiz. Pitch	1.40"	1.63"	1.07"	1.80"	2.75"	2.00"
Label Height	0.50"	0.75"	0.50"	0.75"	1.00"	0.25"
Label Width	1.28"	1.50"	0.94"	1.69"	2.625"	1.50"
No. Across	5	5	7	4	3	4
No. Down	17	12	17	13	10	39
Page Size	8.5" x 11"	8.5" x 11"	8.5" x 11"	8.5" x 11"	8.5" x 11"	8.5" x 11"

See reverse side for more information on Cryo-Babies® and Cryo-Tags® →

LASER CRYO-BABIES[®] AND CRYOTAGS[®]

Labels for Cryogenic Storage*


TOUGH - Chemically inert, polyolefin labels resist most organic solvents and caustic agents. They adhere to most plastics[†], glass, and metals without cracking, peeling, or degrading.

TEMPERATURE-RESISTANT - From -196°C to 100°C, Cryo-Babies and Cryo-Tags withstand boiling water baths, moderate oven temperatures (150°C), autoclaving and extreme freezer temperatures (liquid and vapor phase nitrogen).

SPECIALLY DESIGNED - Label sheets are specially top-coated to readily accept laser printing. Heat-resistant, stay-flat sheets will not jam in laser printers.

IDEALLY SIZED - Will fit a wide variety of cryovials, microtubes, storage boxes and other containers.

COLOR CODING - Each label size is available in White, Blue, Green, Orange, Red and Yellow.

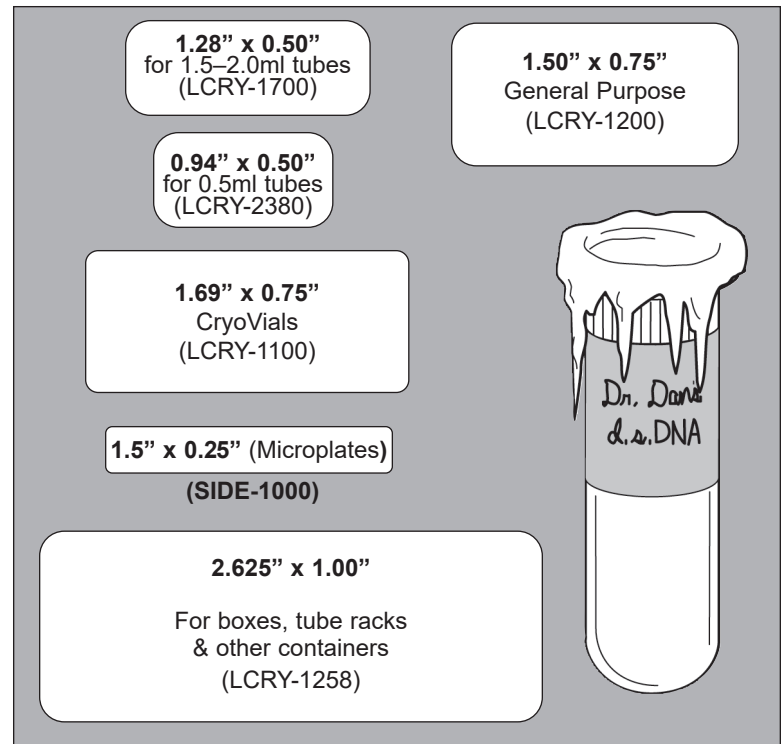

CAT. NO.	COLOR	LABEL SIZE	U OF M
LCRY-1100-(x)	Specify	1.69" x 0.75" / 43mm x 19mm	1,040/pk
LCRY-1200-(x)	Specify	1.50" x 0.75" / 38mm x 19mm	1,200/pk
LCRY-1700-(x)	Specify	1.28" x 0.50" / 33mm x 13mm	1,700/pk
LCRY-2380-(x)	Specify	0.94" x 0.50" / 24mm x 13mm	2,380/pk
SIDE-1000-(x)	Specify	1.50" x 0.25" / 38mm x 6mm	3,120/pk
LCRY-1258-(x)	Specify	2.625" x 1.00" / 67mm x 25mm	600/pk
RNBW-2000	Rainbow*	1.69" x 0.75" / 43mm x 19mm	1,040/pk
RNBW-2100	Rainbow*	1.50" x 0.75" / 38mm x 19mm	1,200/pk
RNBW-2200	Rainbow*	1.28" x 0.50" / 33mm x 13mm	1,700/pk
RNBW-2300	Rainbow*	0.94" x 0.50" / 24mm x 13mm	2,380/pk
RNBW-2400	Rainbow*	1.50" x 0.25" / 38mm x 6mm	3,120/pk
RNBW-2500	Rainbow*	2.625" x 1.00" / 67mm x 25mm	600/pk

To specify color, substitute (x) with: B=Blue, G=Green, O=Orange, R=Red, Y=Yellow
For white, do not include a letter (LCRY-1700 = White)

*Rainbow packs contain 4 sheets each of: Blue, Green, Orange, Red, Yellow

ORDER INFORMATION

Online: www.divbio.com

Phone: 1-800-796-9199
1-781-326-6709

Fax: 1-781-326-5611

Mail: **Diversified Biotech**
65 Commerce Way
Dedham, MA 02026

[†] NOTE: Siliconized surfaces should be avoided. Apply to clean, dry tubes at room temperature. Tubes made of 100% virgin polypropylene are recommended.

*patented